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Deagglomeration and/or sizing machine by millstone type grinder

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Deagglomeration and/or sizing machine by millstone type grinder in elution rate for pharmaceutical use, Supermasscolloider

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不揃いの原薬粒子を効率よく解砕整粒  
Deagglomerate ununiform API particles efficiently.

針状結晶が集積した原薬粒子 (流動不良や二次凝集の原因)  
API crystals have long and thin shapes (Cause of poor flowability or secondary agglomeration)

原薬粒子の季節変動  
Variation of API particles with the seasons

解砕整粒後の原薬  
API particles after deagglomeration and/or sizing

\* API: Active Pharmaceutical Ingredient

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Deagglomeration and/or sizing by grinders

- 大きな粒子は選択的に解砕  
Pulverize larger particle diameters selectively
- 小さい粒子は過粉碎しない  
Small particles are not pulverized excessively
- 静電気帯電が少ない  
Low electrostatic charging

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provides excellence  
PH debut!

Millstone type grinding machine "Supermasscolloider" that has so many experiences with food industry established the technology of deagglomeration and/or sizing for pulverizing, is launched as "Supermasscolloider PH".

This machine will not over pulverize particles depending on different manufactures of API or the seasons, and enable to deagglomeration and/or sizing large particles uniformly.

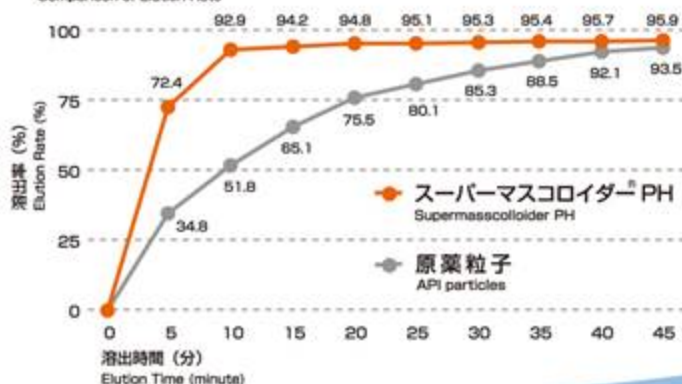
After deagglomeration and/or sizing, the particles will have no secondary agglomeration. The elution rate will dramatically increase and the defective elution rate will dramatically decrease.

Another feature is the cooling effect of air vortex caused by the gap of the grinders, it will cause less heat generation and decomposition of the drug substances.

API particles which have low melting point enable to deagglomeration and/or sizing and has a potential for improving elution rate.

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

Easy to open and close the heavy lid by the spring balancer. It cannot be operated unless assembling all of the parts. Sensors are provided to various places.

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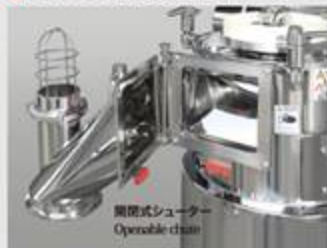Operable chute for ease of maintenance.

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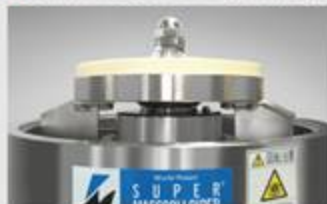Wash to clean the back side of the grinding wheel.

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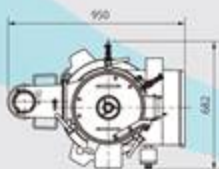
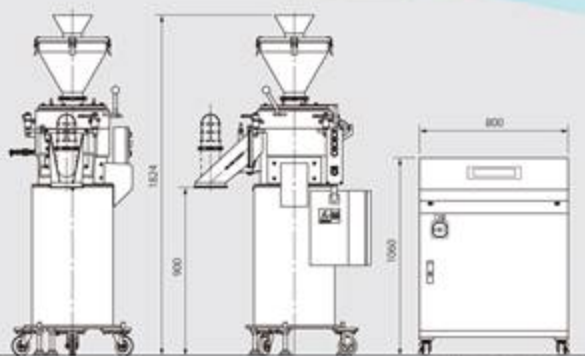
Meister function can be equipped.  
New development "Smart Meister function" can be equipped.  
Meister function provides automatic control of grinder clearance according to preset calibration enabling unattended operation.

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MASUKO SANGYO CO.,LTD.